

TOOLS Prestained Protein Marker 5–245 kDa

Cat. no.: TM-PM05245

Storage: Store at -20°C

Product size: 500 μL

Introduction

TOOLS Pre-Stain Protein marker (large) is ready-to-use and premixed with a loading buffer for direct loading (5 μL) on SDS-PAGE (tris-glycine buffer). This product requires no further dilution or addition of any reducing agent. After the gel is run, the marker yields the following 13 discrete fragments (in kDa): 75 (red), 25 (green), 245, 180, 135, 100, 63, 48, 35, 20, 17, 11, and 5 (blue).

Protocol

1. Load 5 μL for each gel well.
2. Do not boil the marker.

