

TOOLS Prestained Protein Marker large

Cat. no.: TM-PM10315

Storage: Store at -20°C


Product size: 500 μL

Introduction

TOOLS Pre-Stain Protein Marker large is ready-to-use marker, premixed with a loading buffer for direct loading (at 5 μL) on SDS-PAGE (a Tris-glycine buffer). This product requires no further dilution or addition of any reducing agent. After the gel is run, it yields 12 distinct fragments, indicating molecular weights in KDa as follows: 315 (orange), 250 (blue), 180 (blue), 130 (blue), 95 (blue), 72 (orange), 52 (green), 43 (blue), 34 (blue), 26 (blue), 17 (blue), and 10 (green).

Protocol

1. Load 5 μL for each gel.
2. Do not boil the marker.

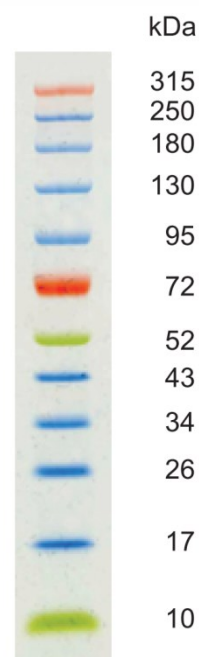

Tris-Glycin
5~20%