

TOOLStripping Buffer

Cat. no.: TW-ST500

Storage: Store at room temperature

Product Size: 500 mL

Introduction

TOOLStripping Buffer is a ready-to-use buffer that safely and effectively removes primary and secondary antibodies from Western blot membranes; simply incubate this buffer with PVDF or a nitrocellulose membrane to allow Western blots to be reprobed for other antibodies.

Protocol

1. Add TOOLStripping Buffer to completely cover the developed membrane and incubate for 3 min at room temperature.
2. Wash the membrane with a large amount of ddH₂O.
3. Wash twice with PBST for 5 min at room temperature.
4. Proceed to reprobe for other antibodies.

Note: TOOLSignal Boosting Buffer (TW-NB250) is recommended for reprobing.

The product is for research use only. It is not for diagnostic or clinical use.

